

ACIDITY ON OIL



Turn on the Minifood.

At the end of the heating, insert the test tube with the reagent in the cell of incubation for 3-5 minut.

At the end of the heating, to do Low Acidity test press **ACID**, press again ACID for High Acidity.



Agitate the test tube by inverting it.

Insert the test tube with reagent in the reading cell and press **ENTER**.



Open the test tube.

Important: Place the cap with the internal part pointing upwards. Do not touch the internal part with the hands.




Agitate the sample several times to make it homogenous.



Taking 2.5 μ l of sample for Low Acidity or 1.5 μ l for High Acidity. Immerse the tip of the pipette into the oil with the piston lowered. **SLOWLY** lift the piston keeping the point immersed in the sample.

Cleaning the tip: let the tip slide onto the paper as indicated in the photograph.

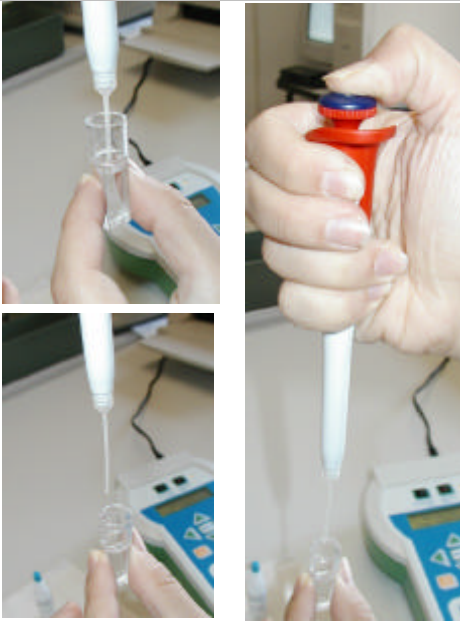

Immerse the tip into the liquid in the test tube and pump the piston up and down 2-3 times by using the thumb.



Plug the test tube and agitate by inverting it.



Insert the test tube in the reading cell and press **ENTER**.



On display you have the result of test.