

PEROXIDES ON OIL



Turn on the Minifood.

At the end of the heating, insert the test tube with the reagent in the cell of incubation for 3-5 minut.

At the end of the heating, to do Peroxides test press **PEROX**.



Agitate the test tube by inverting it.

Insert the test tube with reagent in the reading cell and press **ENTER**.



Open the test tube.

Important: Place the cap with the internal part pointing upwards. Do not touch the internal part with the hands.




Agitate the sample several times to make it homogenous.



Taking 2.5 μ l of sample: immerse the tip of the pipette into the oil with the piston lowered. SLOWLY lift the piston keeping the point immersed in the sample.

Cleaning the tip: let the tip slide onto the paper as indicated in the photograph.

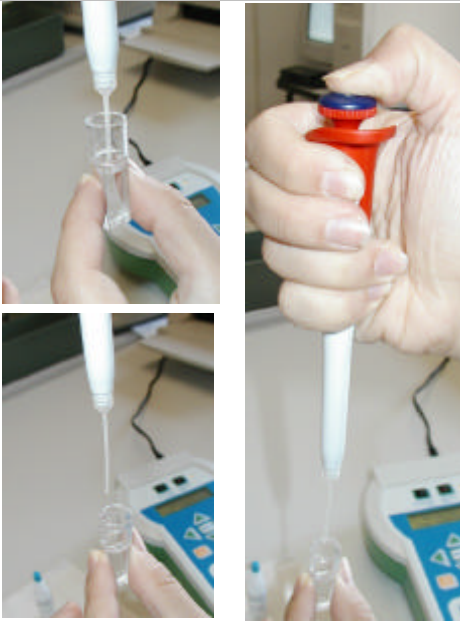

Immerse the tip into the liquid in the test tube and pump the piston up and down 2-3 times by using the thumb.



Plug the test tube and agitate by inverting it.

Open the test tube and insert one drop of R2.

Plug the test tube and agitate by inverting it.



Insert the test tube in the reading cell and press **ENTER**.



After 2 minutes on display you have the result of test.